

OVERVIEW

The Spillchamp 240-Litre Chemical Spill Kit Wheelie Bin is the go-to resource for managing large-scale, aggressive chemical spills in challenging warehouse environments. Built with durability and mobility in mind, it can be swiftly wheeled directly to the spill area, ensuring a fast and effective response.

Inside, the kit is packed with highly absorbent pads and socks specifically designed for aggressive substances like corrosive chemicals, acids, and solvents, offering reliable performance for hazardous cleanups.

Each kit absorbs approximately 240 litres and includes -

- 200 x chemical absorbent pads
- 12 x chemical absorbent sock
- 4 x PVC waste bag and tie
- 1x PPE kit (including goggles and gloves)
- 1x black/yellow barrier tape

PRODUCT FEATURES

- Robust HDPE wheelie bin convenient reusable bin to store and transport the kit
- Large capacity suitable for larger localised chemical spills
- Waste bags and PPE included safe and easy disposal of waste items
- Refillable reusable wheelie bin, refill with spill contents for repeat use
- Suitable for aggressive chemical, acid or solvent spills in labs, warehouses and manufacturing facilities

IRELAND www.pittman.ie • sales@pittman.ie UK www.pittman.uk • sales@pittman.uk

TECHNICAL DATA





SPECIFICATION	CHEMICAL SPILL PADS	CHEMICAL SPILL SOCKS
Material:	Hydrophobic textile	
Color:	Yellow	
Item absorbency:	1 litre	3 - 4 litres
Quantity:	200	12
Length:	500mm	1200mm
Width:	400mm	70mm
Total kit absorbency:	240 litres	
Liquids absorbed:	Acids and aggressive chemicals	
Container type:	HDPE wheelie bin	
Applications	Large capacity suitable for localised aggressive chemical, acid or solvent spills in labs, warehouses and manufacturing facilities	



INSTRUCTIONS FOR USE

- **1. Acquire the spill kit:** Bring the kit to the spill location it is recommended to have the kit close by to places proned to spills
- **2. Cordon spill area:** Prevent access to the area by using barriers or warnings to deter pedestrian entry
- **3. Apply PPE:** Ensure any cleanup staff wear gloves and goggles to avoid injury
- **4. Apply materials to spill:** Apply the pads and socks to the spill and allow to soak as long as necessary.
- **5. Disposal:** Use any included waste bags to dispose of waste material safely and efficiently

